


illumia®Plus2 Integrating Sphere Spectroradiometers

Flexible, modular solutions with centralized data management

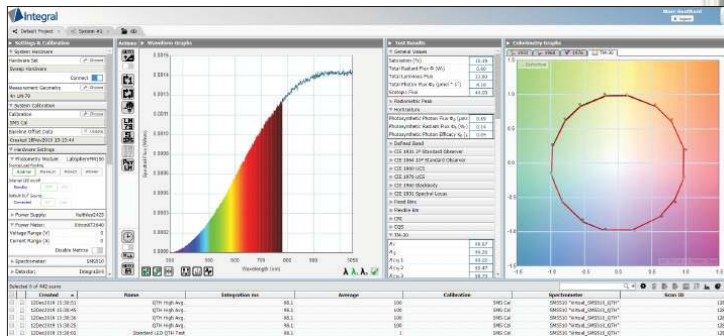

Build a system that meets your needs. Labsphere illumiaPlus2 integrating sphere spectroradiometers offer flexibility to meet your most demanding requirements and grow with your business. A choice of integrating sphere sizes ranging from 25 cm to 3 m, spectrometer options, and application-specific modules help you maximize throughput and productivity while maintaining high absolute accuracy and repeatability.

Integral® Software drives it all with operation from any device, any platform, any location and in any language.

Improve productivity:

- Add-on electronic modules increase functionality and simplify compliance with IES LM-79-19, IES LM-78 and equivalent measurement guidelines
- Automated calibration routines ensure ease-of-use and improved efficiency
- Automated IES LM-79-19 stabilization routine
- Improved integrating sphere design allowing lamp “hot swap” to achieve greater efficiency and shorter measurement time
- Generate reports using Excel templates: data where you want it, how you want it, formatted for language and style

Find the illumiaPlus2 system that best fits your application

illumiaPlus2 800

Measure:

- Packaged LEDs
- Clustered LEDs
- Miniature Lamps
- Entertainment Lighting
- Automotive Lighting
- LED Troffers
- LED Luminaires

Key Features:

- Wide spectral range
- Fast CCD array detector
- Compact size
- Ideal for QC and manufacturing applications
- High sensitivity

illumiaPlus2 2600

Measure:

- Indoor Lighting
- Outdoor Lighting
- Roadway Lighting
- Lamp and Luminaires
- LEDs
- Entertainment Lighting
- Automotive Lighting
- Troffers
- Luminaires
- CFLs
- Fluorescent Lamps
- OLEDs
- Low Power LEDs

Key Features:

- Fast, low noise; TE cooled back – thinned CCD array detector
- Shutter for dark measurements in real time
- Hardware triggering capability
- Exceptional stability at long exposure time
- High dynamic range

illumiaPlus2 3020/3030

Measure:

- Packaged LEDs
- Clustered LEDs
- Miniature Lamps
- Entertainment Lighting
- Automotive Lighting
- LED Troffers
- LED Luminaires
- CFLs
- Fluorescent Lamps
- OLEDs
- Low Power LEDs
- HID Lamps
- LED Devices in High Speed
- Production Environment

Key Features:

- Exceptional measurement sensitivity
- Inbuilt filter wheel with ND filters provides excellent dynamic range
- Low stray light
- High speed
- Easily synchronized with other devices

Every illumiaPlus2 system features these standard products

Programmable DC Power Supplies

Designed to accurately provide DC current to reference lamp, auxiliary lamp, and DCV devices under test.

The current output is selected, set and controlled using Integral Software included with the power supply.

- Programmable regulated DC current
- Programmable regulated DC voltage
- Controlled current ramp up
- Lamp operation timer
- Easy on/off operation
- Front panel or remote control
- Current, voltage readback

ICM-500 Control Module

The illumiaPlus2 Control Module is the routing module that ties Labsphere's powerful Integral Software to the illumiaPlus2 total spectral flux measurement hardware. When the ICM-500 is controlled by Integral, this user friendly, turn key system automatically routes power and metering.

- Main hub for power supplies and power meters
- Routes DC voltage to 2π and 4π reference locations
- Routes power to absorption correction lamp
- Routes DC or AC power to devices under test
- USB inputs




Calibrated Spectral and Luminous Flux Standards

Each standard has been carefully screened, seasoned, and calibrated at our manufacturing facility under the guidelines recommended by the NVLAP accredited ISO 17025 practices for the highest degree of confidence.



Integral® Light Measurement Software

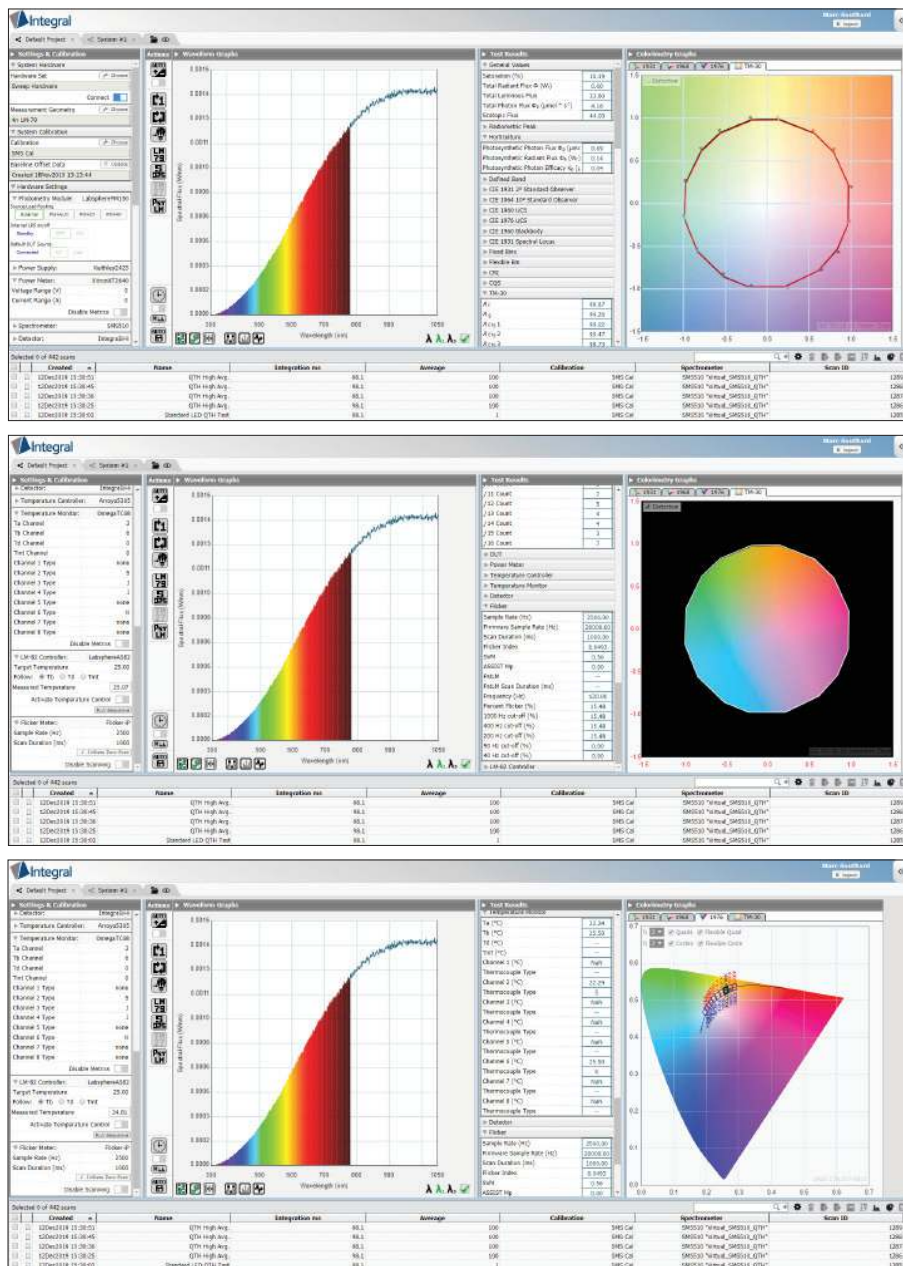

- HTML5-enabled web browser based light measurement software
- Operation from any device, any platform, any location and in any language
- Instantly switch between English, Mandarin Chinese, Japanese, Korean, and French
- Large assortment of test hardware configurations are supported (spectrometer, AC and DC power supplies, temperature controls and monitors)
- Powerful, easy-to-use Application Programming Interface (API) supports LabVIEW, .NET, C, and VBA
- One user can control many test stations and multiple users can access the same test station from anywhere
- Meets LM-79-19 and LM-78 integrating sphere spectrometer recommended measurement methods
- Automated calibration routines
- Built-in report generator with the ability to create custom reports
- All Industry standard color calculations including:
 - x, y, u, v, u', v', CCT, CRI (1-15 and general), CQS, luminous flux (lumens), scotopic lumens, Duv, dominant wavelength, peak wavelength, FWHM, Centroid, Purity, ANSI SSL 2015 binning, TM-30-18 fidelity and gamut data, distortion and vector graphics, and horticulture

Supported Devices		
Spectrometer	Power Supply	Power Meter
CDS1100	Agilent6811B	QINGZH8775C1
CDS1200	Agilent6812B	TETP2201
CDS2400	AgilentE3612A	Xiltron2801
CDS2600	AgilentE3613A	YongguaiWY7210
CDS3000	AgilentE3614A	YongguaiWY73000
CDS3010	AgilentN5751A	YongguaiWY7310
CDS3020	Amptek1501X	
CDS3030	Amptek1501X	
CDS500	Chroma1601	Temperature Controller
CDS610	Chroma1602	Arroyo5300
QEPProGeneric	Chroma1603	Arroyo5305
SMS500	Chroma1604	Arroyo5515
SMS500ULS	Chroma1605	LeindTEC
SMS510	Keithley2400	
	Keithley2410	
	Keithley2420	
	Keithley2425	
	Keithley2430	
	Keithley2440	
	Keithley2450	
	Keithley2460	
	Keithley2470	
	Keithley2480	
	Keithley2490	
	Keithley2500	
	Keithley2510	
	Keithley2520	
	Keithley2530	
	Keithley2540	
	Keithley2550	
	Keithley2560	
	Keithley2570	
	Keithley2580	
	Keithley2590	
	Keithley2600	
	Keithley2610	
	Keithley2620	
	Keithley2630	
	Keithley2640	
	Keithley2650	
	Keithley2660	
	Keithley2670	
	Keithley2680	
	Keithley2690	
	Keithley2700	
	Keithley2710	
	Keithley2720	
	Keithley2730	
	Keithley2740	
	Keithley2750	
	Keithley2760	
	Keithley2770	
	Keithley2780	
	Keithley2790	
	Keithley2800	
	Keithley2810	
	Keithley2820	
	Keithley2830	
	Keithley2840	
	Keithley2850	
	Keithley2860	
	Keithley2870	
	Keithley2880	
	Keithley2890	
	Keithley2900	
	Keithley2910	
	Keithley2920	
	Keithley2930	
	Keithley2940	
	Keithley2950	
	Keithley2960	
	Keithley2970	
	Keithley2980	
	Keithley2990	
	Keithley3000	
	Keithley3010	
	Keithley3020	
	Keithley3030	
	Keithley3040	
	Keithley3050	
	Keithley3060	
	Keithley3070	
	Keithley3080	
	Keithley3090	
	Keithley3100	
	Keithley3110	
	Keithley3120	
	Keithley3130	
	Keithley3140	
	Keithley3150	
	Keithley3160	
	Keithley3170	
	Keithley3180	
	Keithley3190	
	Keithley3200	
	Keithley3210	
	Keithley3220	
	Keithley3230	
	Keithley3240	
	Keithley3250	
	Keithley3260	
	Keithley3270	
	Keithley3280	
	Keithley3290	
	Keithley3300	
	Keithley3310	
	Keithley3320	
	Keithley3330	
	Keithley3340	
	Keithley3350	
	Keithley3360	
	Keithley3370	
	Keithley3380	
	Keithley3390	
	Keithley3400	
	Keithley3410	
	Keithley3420	
	Keithley3430	
	Keithley3440	
	Keithley3450	
	Keithley3460	
	Keithley3470	
	Keithley3480	
	Keithley3490	
	Keithley3500	
	Keithley3510	
	Keithley3520	
	Keithley3530	
	Keithley3540	
	Keithley3550	
	Keithley3560	
	Keithley3570	
	Keithley3580	
	Keithley3590	
	Keithley3600	
	Keithley3610	
	Keithley3620	
	Keithley3630	
	Keithley3640	
	Keithley3650	
	Keithley3660	
	Keithley3670	
	Keithley3680	
	Keithley3690	
	Keithley3700	
	Keithley3710	
	Keithley3720	
	Keithley3730	
	Keithley3740	
	Keithley3750	
	Keithley3760	
	Keithley3770	
	Keithley3780	
	Keithley3790	
	Keithley3800	
	Keithley3810	
	Keithley3820	
	Keithley3830	
	Keithley3840	
	Keithley3850	
	Keithley3860	
	Keithley3870	
	Keithley3880	
	Keithley3890	
	Keithley3900	
	Keithley3910	
	Keithley3920	
	Keithley3930	
	Keithley3940	
	Keithley3950	
	Keithley3960	
	Keithley3970	
	Keithley3980	
	Keithley3990	
	Keithley4000	
	Keithley4010	
	Keithley4020	
	Keithley4030	
	Keithley4040	
	Keithley4050	
	Keithley4060	
	Keithley4070	
	Keithley4080	
	Keithley4090	
	Keithley4100	
	Keithley4110	
	Keithley4120	
	Keithley4130	
	Keithley4140	
	Keithley4150	
	Keithley4160	
	Keithley4170	
	Keithley4180	
	Keithley4190	
	Keithley4200	
	Keithley4210	
	Keithley4220	
	Keithley4230	
	Keithley4240	
	Keithley4250	
	Keithley4260	
	Keithley4270	
	Keithley4280	
	Keithley4290	
	Keithley4300	
	Keithley4310	
	Keithley4320	
	Keithley4330	
	Keithley4340	
	Keithley4350	
	Keithley4360	
	Keithley4370	
	Keithley4380	
	Keithley4390	
	Keithley4400	
	Keithley4410	
	Keithley4420	
	Keithley4430	
	Keithley4440	
	Keithley4450	
	Keithley4460	
	Keithley4470	
	Keithley4480	
	Keithley4490	
	Keithley4500	
	Keithley4510	
	Keithley4520	
	Keithley4530	
	Keithley4540	
	Keithley4550	
	Keithley4560	
	Keithley4570	
	Keithley4580	
	Keithley4590	
	Keithley4600	
	Keithley4610	
	Keithley4620	
	Keithley4630	
	Keithley4640	
	Keithley4650	
	Keithley4660	
	Keithley4670	
	Keithley4680	
	Keithley4690	
	Keithley4700	
	Keithley4710	
	Keithley4720	
	Keithley4730	
	Keithley4740	
	Keithley4750	
	Keithley4760	
	Keithley4770	
	Keithley4780	
	Keithley4790	
	Keithley4800	
	Keithley4810	
	Keithley4820	
	Keithley4830	
	Keithley4840	
	Keithley4850	
	Keithley4860	
	Keithley4870	
	Keithley4880	
	Keithley4890	
	Keithley4900	
	Keithley4910	
	Keithley4920	
	Keithley4930	
	Keithley4940	
	Keithley4950	
	Keithley4960	
	Keithley4970	
	Keithley4980	
	Keithley4990	
	Keithley5000	
	Keithley5010	
	Keithley5020	
	Keithley5030	
	Keithley5040	
	Keithley5050	
	Keithley5060	
	Keithley5070	
	Keithley5080	
	Keithley5090	
	Keithley5100	
	Keithley5110	
	Keithley5120	
	Keithley5130	
	Keithley5140	
	Keithley5150	
	Keithley5160	
	Keithley5170	
	Keithley5180	
	Keithley5190	
	Keithley5200	
	Keithley5210	
	Keithley5220	
	Keithley5230	
	Keithley5240	
	Keithley5250	
	Keithley5260	
	Keithley5270	
	Keithley5280	
	Keithley5290	
	Keithley5300	
	Keithley5310	
	Keithley5320	
	Keithley5330	
	Keithley5340	
	Keithley5350	
	Keithley5360	
	Keithley5370	
	Keithley5380	
	Keithley5390	
	Keithley5400	
	Keithley5410	
	Keithley5420	
	Keithley5430	
	Keithley5440	
	Keithley5450	
	Keithley5460	
	Keithley5470	
	Keithley5480	
	Keithley5490	
	Keithley5500	
	Keithley5510	
	Keithley5520	
	Keithley5530	
	Keithley5540	
	Keithley5550	
	Keithley5560	
	Keithley5570	
	Keithley5580	
	Keithley5590	
	Keithley5600	
	Keithley5610	
	Keithley5620	
	Keithley5630	
	Keithley5640	
	Keithley5650	
	Keithley5660	
	Keithley5670	
	Keithley5680	
	Keithley5690	
	Keithley5700	
	Keithley5710	
	Keithley5720	
	Keithley5730	
	Keithley5740	
	Keithley5750	
	Keithley5760	
	Keithley5770	
	Keithley5780	
	Keithley5790	
	Keithley5800	
	Keithley5810	
	Keithley5820	
	Keithley5830	
	Keithley5840	
	Keithley5850	
	Keithley5860	
	Keithley5870	
	Keithley5880	
	Keithley5890	
	Keithley5900	
	Keithley5910	
	Keithley5920	
	Keithley5930	
	Keithley5940	
	Keithley5950	
	Keithley5960	
	Keithley5970	
	Keithley5980	
	Keithley5990	
	Keithley6000	
	Keithley6010	
	Keithley6020	
	Keithley6030	
	Keithley6040	
	Keithley6050	
	Keithley6060	
	Keithley6070	
	Keithley6080	
	Keithley6090	
	Keithley6100	
	Keithley6110	
	Keithley6120	
	Keithley6130	
	Keithley6140	
	Keithley6150	
	Keithley6160	
	Keithley6170	
	Keithley6180	
	Keithley6190	
	Keithley6200	
	Keithley6210	
	Keithley6220	
	Keithley6230	
	Keithley6240	
	Keithley6250	
	Keithley6260	
	Keithley6270	
	Keithley6280	
	Keithley6290	
	Keithley6300	
	Keithley6310	
	Keithley6320	
	Keithley6330	
	Keithley6340	
	Keithley6350	
	Keithley6360	
	Keithley6370	
	Keithley6380	
	Keithley6390	
	Keithley6400	
	Keithley6410	
	Keithley6420	
	Keithley6430	
	Keithley6440	
	Keithley6450	
	Keithley6460	
	Keithley6470	
	Keithley6480	
	Keithley6490	
	Keithley6500	
	Keithley6510	
	Keithley6520	
	Keithley6530	
	Keithley6540	
	Keithley6550	
	Keithley6560	
	Keithley6570	
	Keithley6580	
	Keithley6590	
	Keithley6600	
	Keithley6610	
	Keithley6620	
	Keithley6630	
	Keithley6640	
	Keithley6650	
	Keithley6660	
	Keithley6670	
	Keithley6680	
	Keithley6690	
	Keithley6700	
	Keithley6710	
	Keithley6720	
	Keithley6730	
	Keithley6740	
	Keithley6750	
	Keithley6760	
	Keithley6770	
	Keithley6780	
	Keithley6790	
	Keithley6800	
	Keithley6	

illumiaPlus2 - 800 Ordering Information

System:	illumiaPlus2 with 25 cm sphere	illumiaPlus2 with 50 cm sphere	illumiaPlus2 with 100 cm sphere	illumiaPlus2 with 165 cm sphere	illumiaPlus2 with 195 cm sphere
illumiaPlus2 800					
2 π version:	—	AA-40022-050	AA-40022-100	AA-40022-165	AA-40022-195
4 π version:	AA-40024-025	AA-40024-050	AA-40024-100	AA-40024-165	AA-40024-195
Above Systems Include:					
Sphere:	LMS-025	LMS-050	LMS-100	LMS-165	LMS-195
Control Module:	ICM-500	ICM-500	ICM-500	ICM-500	ICM-500
DC Power Supply:	LPS-175-27	LPS-175-27	LPS-525-42	LPS-525-42	LPS-525-42
Spectrally-Calibrated Lamp: (2 π and 4 π are not interchangeable)					
2 π	2PI-1-INT-650	2PI-1-INT-650	2PI-1-INT-1400	2PI-1-INT-1400	2PI-1-INT-1400
4 π	SCL-650	SCL-650	SCL-1400	SCL-1400	SCL-1400
Aux Lamp:	AUX-650	AUX-650	AUX-1400	AUX-1400	AUX-1400
Lamp Socket Assembly:					
	2 π /4 π	2 π /4 π	2 π /4 π	2 π /4 π	2 π /4 π
Software:	Integral	Integral	Integral	Integral	Integral

All systems include Retouch 6080 and Tool Kit

Typical Performance Specifications (lumens)

Spectral Flux Measurements: 350 nm - 1000 nm

Exposure Time Range: 40 μ s – 5 sec (Actual exposure time depends on sphere size and source type)

	min	max	min	max	min	max	min	max	min	max
Tungsten Filament:	0.039	1880	0.154	7500	0.616	16300	1.39	41000	2.46	57000
Cool White LED:	0.001	1770	0.003	7100	0.013	14500	0.030	37000	0.054	52000
Warm White LED:	0.001	1100	0.005	4500	0.020	13800	0.045	34000	0.080	47000
Blue LED:	0.025	77	0.099	300	0.398	1200	0.895	2700	1.59	3800
Red LED:	0.009	195	0.037	800	0.148	1100	0.334	3500	0.594	5000
Upper Range:	Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C	

illumiaPlus2 - 2600 Ordering Information

System:	illumiaPlus2 with 25 cm sphere	illumiaPlus2 with 50 cm sphere	illumiaPlus2 with 100 cm sphere	illumiaPlus2 with 165 cm sphere	illumiaPlus2 with 195 cm sphere
illumiaPlus2 2600					
2π version:	—	AA-40072-050	AA-40072-100	AA-40072-165	AA-40072-195
4π version:	AA-40074-025	AA-40074-050	AA-40074-100	AA-40074-165	AA-40074-195
Above Systems Include:					
Sphere:	LMS-025	LMS-050	LMS-100	LMS-165	LMS-195
Control Module:	ICM-500	ICM-500	ICM-500	ICM-500	ICM-500
DC Power Supply:	LPS-175-27	LPS-175-27	LPS-525-42	LPS-525-42	LPS-525-42
Spectrally-Calibrated Lamp: (2π and 4π are not interchangeable)					
2π	2PI-1-INT-650	2PI-1-INT-650	2PI-1-INT-1400	2PI-1-INT-1400	2PI-1-INT-1400
4π	SCL-650	SCL-650	SCL-1400	SCL-1400	SCL-1400
Aux Lamp:	AUX-650	AUX-650	AUX-1400	AUX-1400	AUX-1400
Lamp Socket Assembly:					
	2π/4π	2π/4π	2π/4π	2π/4π	2π/4π
Software:	Integral	Integral	Integral	Integral	Integral

All systems include Retouch 6080 and Tool Kit

Typical Performance Specifications (lumens)

Spectral Flux Measurements: 350 nm - 1050 nm

Exposure Time Range: 8 ms - 900 sec (Actual exposure time depends on sphere size and source type)

	min	max	min	max	min	max	min	max	min	max
Tungsten Filament:	0.007	3400	0.031	13800	0.12	22000	0.33	89000	0.47	124000
Cool White LED:	0.003	1900	0.012	7600	0.04	30000	0.13	114000	0.18	160000
Warm White LED:	0.002	1300	0.009	5300	0.03	25000	0.09	98500	0.14	138000
Blue LED:	0.001	140	0.002	600	0.01	1800	0.02	6000	0.04	8300
Red LED:	0.002	330	0.007	1300	0.03	1900	0.07	7600	0.11	11000
Upper Range:	Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C	

illumiaPlus2 - 3020/3030 Ordering Information

System:	illumiaPlus2 with 50 cm sphere	illumiaPlus2 with 100 cm sphere	illumiaPlus2 with 165 cm sphere	illumiaPlus2 with 195 cm sphere
illumiaPlus2 3020:				
2 π version:	AA-40082-050	AA-40082-100	AA-40082-165	AA-40082-195
4 π version:	AA-40084-050	AA-40084-100	AA-40084-165	AA-40084-195
illumiaPlus2 3030:				
2 π version:	AA-40092-050	AA-40092-100	AA-40092-165	AA-40092-195
4 π version:	AA-40094-050	AA-40094-100	AA-40094-165	AA-40094-195
Above Systems Include:				
Sphere:	LMS-050	LMS-100	LMS-165	LMS-195
Control Module:	ICM-500	ICM-500	ICM-500	ICM-500
DC Power Supply:	LPS-175-27	LPS-175-27	LPS-175-27	LPS-175-27
Spectrally-Calibrated Lamp: (2 π and 4 π are not interchangeable)				
2 π :	2PI-1-INT-040	2PI-1-INT-040	2PI-1-INT-650	2PI-1-INT-650
4 π :	SCL-040	SCL-040	SCL-650	SCL-650
Aux Lamp:	AUX-040	AUX-040	AUX-650	AUX-650
Lamp Socket Assembly:	2 π /4 π	2 π /4 π	2 π /4 π	2 π /4 π
Software:	Integral	Integral	Integral	Integral

All systems include Retouch 6080 and Tool Kit

Typical Performance Specifications (lumens)

Spectral Flux Measurements: 350 nm - 830 nm (3020 systems) 350 nm - 1050 nm (3030 systems)

Exposure Time Range: 5 ms - 20 sec (Actual exposure time depends on sphere size and source type)

	min	max	min	max	min	max	min	max
Tungsten Filament:	0.025	19000	0.100	76000	0.26	210000	0.37	291000
Cool White LED:	0.007	20800	0.013	83000	0.03	227000	0.05	317000
Warm White LED:	0.004	20000	0.009	79000	0.02	215000	0.03	300000
Blue LED:	0.010	1100	0.018	4541	0.05	12300	0.07	17000
Red LED:	0.008	1600	0.015	6245	0.04	17000	0.05	24000
Upper Range:	Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C	

illumiaPlus2 - 3 Meter Light Measurement Systems Ordering Information

System:	illumiaPlus2 600-300	illumiaPlus2 610-300	illumiaPlus2 2600-300	illumiaPlus2 3020-300	illumiaPlus2 3030-300
4π version:	AA-40024-300	AA-40034-300	AA-40074-300	AA-40084-300	AA-40094-300
Sphere:	LMS-300	LMS-300	LMS-300	LMS-300	LMS-300
Control Module:	ICM-500	ICM-500	ICM-500	ICM-500	ICM-500
Power Supply:	LPS-525-42	LPS-525-42	LPS-525-42	LPS-525-42	LPS-525-42
Spectrometer:	CDS-800	CDS-800	CDS-2600	CDS-3020	CDS-3030
Spectrally-Calibrated Lamp:					
4 π :	SCL-1400	SCL-1400	SCL-1400	SCL-1400	SCL-1400
Aux Lamp:	AUX-1400	AUX-1400	AUX-1400	AUX-1400	AUX-1400
Lamp Socket Assembly:	4 π	4 π	4 π	4 π	4 π
Software:	Integral LM	Integral LM	Integral LM	Integral LM	Integral LM

All systems include Retouch 6080 and Tool Kit

Sphere Weight:	1020 lbs (462 kg)
Sphere Max Height:	140 inches (3.6 cm)
Sphere Foot Print Open/Closed:	150 in W x 201 L

Typical Performance Specifications (lumens)

	illumiaPlus2 800-300		illumiaPlus2 2600-300		illumiaPlus2 3020-300 & 3030-300	
	min	max	min	max	min	max
Tungsten Filament:	5.54	128250	1.06	279000	0.83	654750
Cool White LED:	0.12	117000	0.41	360000	0.11	713250
Warm White LED:	0.18	105750	0.32	310500	0.07	675000
Blue LED:	3.58	8550	0.09	18675	0.16	38250
Red LED:	1.34	11250	0.25	24750	0.11	54000
Upper Range:	Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C		Ambient temp can not exceed 100°C	

Integrating Sphere Specifications

Model Number:	LMS-025	LMS-050	LMS-100	LMS-165	LMS-195
Order Number:	AA-30000-025	AA-30000-050	AA-30000-100	AA-30000-165	AA-30000-195
Sphere Size:	25 cm	50 cm	100 cm	165 cm	195 cm
Required Footprint:					
closed	45.9 cm x 30.2 cm x 39.1 cm	74.1 cm x 50.4 cm x 74.0 cm	128.0 cm x 98.1 cm x 166.7 cm	185 cm x 165 cm x 200 cm	213.2 cm x 187 cm x 226 cm
open	64.2 cm x 46.8 cm x 39.1 cm	103.7 cm x 89.9 cm x 74.0 cm	174.4 cm x 165.9 cm x 166.7 cm	269.3 cm x 271.8 cm x 200 cm	310 cm x 307.4 cm x 226 cm
Sphere Weight:	4 Kg	20 Kg	92 Kg	300 Kg	347 Kg
Sphere Color:	Labsphere Blue	Labsphere Blue	Labsphere Blue	Labsphere Blue	Labsphere Blue
Sphere Open Style:	Clam Shell	Clam Shell	Clam Shell	Clam Shell	Clam Shell
Sphere Assembly:	Spun Aluminum	Spun Aluminum	Spun Aluminum	Spun Aluminum	Spun Aluminum
Frame Style:	Extruded Aluminum	Extruded Aluminum	Extruded Aluminum	Extruded Aluminum	Extruded Aluminum
Sphere Coating:	Spectrafect	Spectrafect	Spectrafect	Spectrafect	Spectrafect
Spectrafect Coating Reflectance:	> 97% (nominal)	> 97% (nominal)	> 97% (nominal)	> 97% (nominal)	> 97% (nominal)
Total Number of Ports:	6	6	6	6	6
SMA Adapter:	Included	Included	Included	Included	Included
Cosine Corrector:	Included	Included	Included	Included	Included
External DUT 2π Port Size:	5.08 cm	15 cm	33 cm	53 cm	63 cm
Detector Port Dimension:	1.25 cm	1.25 cm	1.25 cm	1.25 cm	1.25 cm
Detector Port Quantity:	1	2	2	2	2
Port Frame Reducer:	n/a	15 - 2.5 cm 15 cm - 2.5 cm	33 cm - 15 cm; 15 cm - 2.5 cm	53 cm - 15 cm; 15 cm - 2.5 cm	63 cm - 15 cm;
Temperature Probe Port:	2.5 cm	2.5 cm	2.5 cm	2.5 cm	2.5 cm
Max Recommended Lamp Size: (LM-79) 4π geometry	2.5 cm	<7 cm dia, 33 cm long	<14 cm dia, 67 cm long	<23 cm dia, 110 cm long	<27 cm dia, 130 cm long
Max Port or DUT Area: (2π)	5.08 cm dia.; 20 cm ²	15 cm dia.; 71 cm ²	33 cm dia.; 337 cm ²	53 cm dia.; 879 cm ²	63 cm dia.; 1246 cm ²
Max Recommended Linear DUT Dimension: (2/3 sphere diameter)	15 cm	33 cm	66 cm	110 cm	130 cm
Max Recommended Internal DUT Surface Area: (2% Rule, LM-79)	39 cm	157 cm	628 cm	1711 cm	2389 cm
Maximum Sphere Coating Temp:	100°C	100°C	100°C	100°C	100°C

Upgrade Modules Ordering Information

Model Number: IL-AC1

Order Number: AS-40000-002

Includes:

- Chroma 61603 Programmable Instrument Grade AC Power Source
- Cabling for ICM-500 connections

Model Number: IL-PM1

Order Number: AS-40000-001

Includes:

- XITRON 2640 Precision Multi-Channel Power Analyzer
- Cabling for ICM-500 and AC power source connections

Optional Accessories Ordering Information

illumiaPlus2 - 800 and 2600

	25 cm sphere	50 cm sphere	100 cm sphere	165 cm sphere	195 cm sphere
4 π Kit (To perform 4 π measurement geometry with 2 π systems. Includes lamp post, baffle, junction box, SCL calibrated in the 4 π geometry)					
Order Number:	Included	AA-30001-050	AA-30001-100	AA-30001-165	AA-30001-195

4 π Kit (To perform 4 π measurement geometry with 2 π systems. Includes lamp post, baffle, junction box, does not include calibration lamp)					
Order Number:	Included	AA-30006-050	AA-30006-100	AA-30006-165	AA-30006-195

Ambient Temperature Probe and Monitor

Model Number:	TPM-100TC-08	TPM-200TC-08	TPM-400TC-08	TPM-400TC-08	TPM-400TC-08
Order Number:	AS-03003-100	AS-03003-200	AS-03003-400	AS-03003-400	AS-03003-400

Forward Flux Standard

Model Number:	FFS-100-400	FFS-100-400	FFS-100-1000	FFS-100-1000	FFS-100-1000
Order Number:	AS-02768-100	AS-02768-100	AS-02768-200	AS-02768-200	AS-02768-200

illumiaPlus2 - 3020/3030

	50 cm sphere	100 cm sphere	165 cm sphere	195 cm sphere
4 π Kit (To perform 4 π measurement geometry with 2 π systems. Includes lamp post, baffle, junction box, SCL calibrated in the 4 π geometry)				
Order Number:	AA-30005-050	AA-30005-100	AA-30005-165	AA-30005-195

4 π Kit (To perform 4 π measurement geometry with 2 π systems. Includes lamp post, baffle, junction box, does not include calibration lamp)				
Order Number:	AA-30006-050	AA-30006-100	AA-30006-165	AA-30006-195

Ambient Temperature Probe and Monitor

Model Number:	TPM-200TC-08	TPM-400TC-08	TPM-400TC-08	TPM-400TC-08
Order Number:	AS-03003-200	AS-03003-400	AS-03003-400	AS-03003-400

Forward Flux Standard

Model Number:	FFS-100-400	FFS-100-1000	FFS-100-1000	FFS-100-1000
Order Number:	AS-02768-100	AS-02768-200	AS-02768-200	AS-02768-200

illumiaPlus to illumiaPlus2 Upgrade Kit Ordering Information

Model Number: ICM-500-175

Order Number: AS-40000-175

includes: ICM-500, LPS-175 27 DC Power Supply, jumper cable and documentation for systems using 2PI-INT-050, 2PI-INT-650, SCL-040, SCL-650, AUX-040, AUX-650 and FFS-100-400 lamps

Model Number: ICM-500-350

Order Number: AS-40000-350

includes: ICM-500, LPS-350 28 DC Power Supply, jumper cable and documentation for systems using AUX-75, FFS-100-1000, and AUX-100 lamps

Model Number: ICM-500-525

Order Number: AS-40000-525

includes: ICM-500, LPS-525 42 DC Power Supply, jumper cable and documentation for systems using 2PI-INT-1400, AUX-1400, ISC-1400, and SCL-1400 lamps

Model Number: Integral LM-User ASM

Order Number: AS-81021-000

Integral Major Module Software Upgrade to existing illumia and/or Integral installation. Single user, single Integral License and 1 year support and maintenance

